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Product Manual

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Buffered Listeria Enrichment Broth Supplement

Product No.	Product Type	Specification
SR0970	Reagent A: Freeze dried reagent	2*5pieces/box
	Reagent B: Freeze dried reagent	

Intended Use

Used for selective enrichment culture for *Listeria monocytogenes*.

Formula

Product Code	Ingredients	/piece
SR0970A	Nalidixic acid	4mg
	Amphotericin B	1mg
SR0970B	Acridine hydrochloride	1.5mg

Directions

- Opening method of vial: Before opening the vial, the surface of the vial was disinfected with 75 % alcohol cotton. Under sterile conditions, the aluminum cap was opened according to the direction of the arrow on the aluminum cap, and the aluminum cap was torn open the vial rubber stopper.
- Usage of reagent: Each vial was added with 2mL sterile deionized water to completely dissolve the reagent. Add 1 vial each of Listeria Selective Enrichment Supplement (SR0970A and SR0970B) to every 500 mL of the basal medium (HCM190) to prepare Buffered Listeria Enrichment Broth.

Storage and Shelf Life

2~8°C, shelf life on the packaging box

Precautions

- When weighing the dehydrated medium, please wear masks to avoid causing respiratory system discomfort
- Keep container tightly closed after using to prevent clumping.

Waste Disposal

Microbiological contamination was disposed by autoclaving at 121°C for 30 minutes.

Revision

On Dec.1, 2025

GUANGDONG HUANKAI MICROBIAL SCI.&TECH.CO.,LTD.

Address:No. 788 Shenzhou Road, Science City, Huangpu District, Guangzhou, Guangdong, China

Email:support@hkmbio.com

Website:https://www.huankaigroup.com